



220A NAALI



MP-SPR NAVI™ 220A NAALI



MP-SPR Navi™ 220A NAALI provides premium quality data with PureKinetics™ and it also allows measurements of conformation changes.

www.bionavis.com



C3 PROZESS- UND
ANALYSENTECHNIK

MP-SPR NAVI™ 220A NAALI SPECIFICATIONS

Measurement principle	Real-time and label-free Multi-Parametric Surface Plasmon Resonance (MP-SPR) is based on a true goniometric SPR arrangement with a rotating detector. Exceptionally wide angular range 40-78° is measured, real angular resolution 0.001°.
Liquid handling	Automated 96/384-well plate liquid handling for unattended runs . Controlled buffer flow conditions with precise syringe pumps and integrated degasser. 2 separate flow channels . Software switch to select serial or parallel injection mode. Flow rate range from 1 µl/min up to 1000 µl/min.
Sample consumption	Partial loop injections enabling minimized sample consumption. Minimum sample volume 70 µl. Minimum injected volume 50 µl.
Wavelength of light	Standard 670 nm and 785 nm lasers in both flow channels. With additional options, each flow channel can be equipped with 3 lasers (670, 785 and 980nm) or 4 lasers (670, 785, 850 and 980nm). Additional lasers are utilized to detect conformation changes and allow layer characterization.
Refractive index range	1.00-1.40 (measurement bulk environment) which can be extended with additional wavelengths. Layers that MP-SPR can measure may have much higher RI, such as diamond like carbon (2.7) and inorganic crystals. MP-SPR determines also complex refractive index.
Media	One scan encompasses both environments: gas and liquid . Measure not only in water based liquids but also in some organic liquids , such as ethanol and acetonitrile. PureKinetics™ enables liquid composition correction between running buffer and the sample (bulk effect).
Mode of operation	Angular Scanning mode, or "MP-SPR mode": scanning across a range of angles providing complete SPR curve and multiple parameters. Several types of sensograms can be extracted from the recorded SPR curves, such as PureKinetics™. Sampling rate depends selected angular range and resolution, typically less than 2 seconds. Fixed Angle mode, or "traditional SPR mode": measurement from single angle, providing time – intensity sensogram. Mode for fast kinetic studies - sampling rate from 1 ms.
Measurement range	Kinetics: $k_a = 10^3 - 10^8$ 1/(M*s), $k_d = 10^{-7} - 0.1$ 1/s, $K_D = 10^{-3} - 10^{-12}$ M Thickness from Ångströms to micrometers (true range depends on refractive index of the material).
Sensitivity	Smallest detected molecule: In gas: Hydrogen (2 Da); in liquid: small molecules <100 Daltons
Noise	Short-term noise 0.3 µRIU, Baseline drift (long term) < 1µRIU/min.
Temperature	Measurement temperature range from 15 – 45 °C (7° below to 20°C above ambient). Well plate temperature range from 4-40 °C.
Prism	Prism with elastomer enables quick sensor exchange and enables further analysis of sample with other methods.
Flow cells	Standard: 2 channel PDMS flow cell. Flow cell volume is 1 µl. Optional flow cells: SPR321-EC: Electrochemical flow cell, SPR305-MS: High chemical resistant PEEK flow cell, SPR310-GS: Gas flow cell, SPR302-LS: for fast kinetic studies, SPR303-C: Cell flow cell, SPR323-F3: Fluorescence flow cell or request custom made flow cell.
Sensors/Substrates	Wide range of sensor surfaces is available, such as metals (Au, Ag, Cu, Pt etc.), other inorganics (SiO ₂ , Al ₂ O ₃ , TiO ₂ etc.) or functionalized surfaces (CMD, Ni ²⁺ etc.). The sensor holder allows for easy <i>ex-situ</i> depositions on sensors (e.g. spin coating). If you cannot find what you are looking for, request a custom made surface!
Software	Unlimited MP-SPR Navi™ Control and DataViewer software. TraceDrawer™ software for kinetic analysis included. Optional: LayerSolver™ fitting tool for layer characterization.
Maintenance	No mandatory service contract required. Fluidic parts are easily exchanged. Annual maintenance contract (AMC) recommended.
Computer requirements	Win 10 or Win 11, 1 x USB 2.0, 4GB RAM, 10GB hard disk space (1GB for installation + space for measured data)
Dimensions & Weight	W 62 x H 41 x D 47 cm (24" x 16" x 18.5"), 50 kg (110 lbs)
Power requirements	100-240V, 50/60Hz, max. 100W

Specifications are subject to change without prior notice.

Information in this catalogue is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions.



**C3 PROZESS- UND
ANALYSENTECHNIK**

Peter-Henlein-Str. 20
D-85540 Haar b. München
Telefon 089/45 60 06 70
Telefax 089/45 60 06 80
info@c3-analysentechnik.de
www.c3-analysentechnik.de



MP-SPR Navi™ 220A NAALI