



420A ILVES



MP-SPR NAVI™ 420A ILVES



4-channel MP-SPR Navi™ 420A ILVES for molecular interaction studies is the fastest member of the MP-SPR family! Kinetic Titration function makes it an ideal solution for challenging-to-regenerate surfaces, or in case you just want to obtain your results faster!

www.bionavis.com



C3 PROZESS- UND
ANALYSENTECHNIK

MP-SPR NAVI™ 420A ILVES SPECIFICATIONS

Measurement principle	Real-time and label-free Multi-Parametric Surface Plasmon Resonance (MP-SPR) is based on a true goniometric SPR arrangement with a rotating detector. Exceptionally wide angular range 40-78° is measured, real angular resolution 0.001°.
Liquid handling	Automated 96/384-well plate liquid handling for unattended runs . 4 flow channels . Configure channels one-by-one (individually) or in series. Dual loop injection mode enabling e.g. fast Kinetic Titration experiments, and experiments in multiple buffers. Controlled buffer flow conditions with precise syringe pumps and integrated degasser . Flow rate range from 1 µl/min up to 1000 µl/min.
Sample consumption	Partial loop injections enabling minimized sample consumption. Minimum sample volume 70 µl. Minimum injected volume 50 µl.
Wavelength of light	Standard 670 nm and 785nm lasers in all four flow channels. Other wavelengths available on request. Additional wavelengths enable measurement of thickness and refractive index simultaneously.
Refractive index range	1.00-1.40 (measurement bulk environment) which can be extended with additional wavelengths. Layers that MP-SPR can measure may have much higher RI such as diamond-like carbon (2.7) and inorganic crystals. MP-SPR determines also complex refractive index.
Media	One scan encompasses both environments: gas and liquid . Measure not only in water based liquids but also in some organic liquids , such as ethanol and acetonitrile. PureKinetics™ enables liquid composition correction between running buffer and the sample (bulk effect).
Mode of operation	Angular Scanning mode, or "MP-SPR mode": scanning across a range of angles providing complete SPR curve and multiple parameters. Several types of sensograms can be extracted from the recorded SPR curves, such as PureKinetics™. Sampling rate depends selected angular range and resolution, typically less than 2 seconds. Fixed Angle mode, or "traditional SPR mode": measurement from single angle, providing time – intensity sensogram. Mode for fast kinetic studies - sampling rate from 1 ms.
Measurement range	Kinetics: $k_a = 10^3 - 10^8 \text{ 1/(M*s)}$, $k_d = 10^{-7} - 0.1 \text{ 1/s}$, $K_D = 10^{-3} - 10^{-12} \text{ M}$ Thickness from Ångströms to micrometers (true range depends on refractive index of the material).
Sensitivity	Smallest detected molecule: In gas: Hydrogen (2 Da); in liquid: small molecules <100 Daltons
Noise	Short-term noise 0.3 µRIU, Baseline drift (long term) < 1µRIU/min.
Temperature	Measurement temperature range from 15 – 45 °C (7° below to 20°C above ambient). Well plate temperature range from 4-40 °C.
Prism	Prism with elastomer enables quick sensor exchange and enables further analysis of sample with other methods.
Flow cells	Standard: 4 channel PDMS flow cell. Flow cell volume: 1 µl. Optional flow cells: SPR321-EC: Electrochemical flow cell (one chamber), SPR343-C: Cell flow cell or request custom made flow cell.
Sensors/Substrates	Wide range of sensor surfaces is available, such as metals (Au, Ag, Cu, Pt etc.), other inorganics (SiO ₂ , Al ₂ O ₃ , TiO ₂ etc.) or functionalized surfaces (CMD, Ni ²⁺ etc.). The sensor holder allows for easy <i>ex-situ</i> depositions on sensors (e.g. spin coating). If you cannot find what you are looking for, request a custom made surface!
Software	Unlimited MP-SPR Navi™ Control and DataViewer software. TraceDrawer™ software for kinetic analysis and LayerSolver™ fitting tool for layer characterization included.
Maintenance	No mandatory service contract required. Fluidic parts are easily exchanged. Annual maintenance contract (AMC) recommended.
Computer requirements	Win 10 or Win 11, 1 x USB 2.0, 4GB RAM, 10GB hard disk space (1GB for installation + space for measured data)
Dimensions & Weight	W 62 x H 41 x D 47 cm (24" x 16" x 18.5"), 50 kg (110 lbs)
Power requirements	100-240V, 50/60Hz, max. 100W

Specifications are subject to change without prior notice.

Information in this catalogue is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions.



C3 PROZESS- UND
ANALYSENTECHNIK

Peter-Henlein-Str. 20
D-85540 Haar b. München
Telefon 089/45 60 06 70
Telefax 089/45 60 06 80
info@c3-analysentechnik.de
www.c3-analysentechnik.de



MP-SPR Navi™ 420A ILVES